



**For research use only, not for use in diagnostic procedures.**  
Not available for clinical diagnostic use in the United States.

Purification steps
12. Place the tube in the microcentrifuge in a known orientation. Centrifuge at room temperature for 2 minutes at 13,000 rpm (15,000 × <i>g</i> ).
13. Carefully remove the supernatant with a pipette tip and discard it. Take care to avoid disturbing the DNA pellet.
14. Carefully add 250 µL of 70% ethanol. Let stand at room temperature for 1 minute. Completely remove the ethanol without disturbing the pellet.
15. Add 100 µL of TE solution to dissolve the DNA pellet. Vortex for at least 5 seconds.

**Technical support is available Monday to Friday (9h00 to 17h00 ET):**

- Toll-free (North America): 1.866.813.6354, option 6
- All other countries: +1.613.723.5757, option 6
- Email: [support@dnagenotek.com](mailto:support@dnagenotek.com)

Some DNA Genotek products may not be available in all geographic regions.

\*OMNIgene and prepIT are registered trademarks of DNA Genotek Inc. All other brands and names contained herein are the property of their respective owners. All DNA Genotek protocols, white papers and application notes, are available in the support section of our website at [www.dnagenotek.com](http://www.dnagenotek.com).

