


REF OGR-500

**CEF OGR-600** 

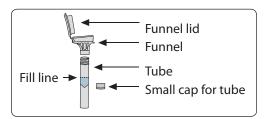

### **Collection precautions:**

Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.

Do NOT remove the plastic film from the funnel lid.

This product is designed for the collection of human DNA from saliva samples.

Contents: Kit contains stabilizing liquid.

**Warnings and precautions:** Wash with water if stabilizing liquid comes in contact with eyes or skin. Do not ingest. See MSDS at www.dnagenotek.com.

Small cap, may pose a choking hazard.

Storage: 15°C 1/30°C

# Summary and explanation of the kit:

Oragene DISCOVER is a self-collection kit that provides the materials and instructions for collecting and stabilizing saliva specimens.

## **Label legend:**

Consult package insert
Collect saliva by (Use by)
REF Catalog number

15°C/t 30°C Storage instructions

Manufacturer

Lot number

SN Serial number

# **USER INSTRUCTIONS**

## Read all instructions prior to collection

### **Procedure:**


Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 5.




Spit into funnel until the amount of liquid saliva (not bubbles) reaches the fill line shown in picture #1.



Hold the tube upright with one hand. Close the funnel lid with the other hand (as shown) by firmly pushing the lid until you hear a loud click. The liquid in the lid will be released into the tube to mix with the saliva. Make sure that the lid is closed tightly.

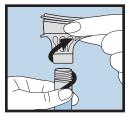

Hold the tube upright. Unscrew the funnel from the tube.

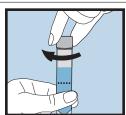

Use the small cap to close the tube tightly.



Shake the capped tube for 5 seconds.
Discard or recycle the funnel.

#### For Research Use Only Not for use in diagnostic procedures

