DNAgenotek

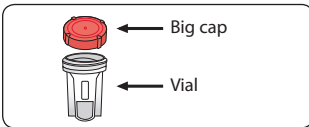
Made in Canada
DNA Genotek Inc.
3000 - 500 Palladium Drive
Ottawa, ON, Canada K2V 1C2

Tel.: +1.613.723.5757
info@dnagenotek.com
www.dnagenotek.com

*Superior samples
Proven performance*

ENGLISH User instructions

Read all instructions prior to collection.

This product is designed for the collection of human RNA from saliva samples.

Contents: Kit contains stabilizing liquid.

Warnings and precautions: Wash with water if stabilizing liquid comes in contact with eyes or skin. Do not ingest. See MSDS at www.dnagenotek.com.

Storage: 15°C / 30°C

Summary and explanation of the kit:

Oragene-RNA is a self-collection kit that provides the materials and instructions for collecting and stabilizing saliva specimens.

Collection precautions:

Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.

Do NOT remove the plastic film from the cap.

Procedure:

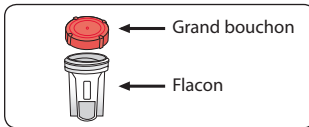
Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 4.

- As saliva gathers in your mouth, begin giving your sample.
- Spit until the amount of liquid saliva (not bubbles) reaches the level shown in picture #2.
- Hold the vial upright. Carefully cap the vial and tighten firmly as shown in picture #3. The stabilizing liquid will be released from the cap.
- Shake the capped vial vigorously for 10 seconds to mix saliva with the liquid from the cap. Discard or recycle outer packaging.

Label legend:

- Consult package insert
- Collect saliva by (Use by)
- Catalog number
- Caution, consult instructions for use
- Storage instructions
- Manufacturer
- Lot number

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all rights reserved.
© 2016 DNA Genotek Inc.,
a subsidiary of OraSure Technologies, Inc.,