

For the collection and stabilization of human and microbial DNA and RNA

Contents: Kit contains stabilizing liquid.

Warnings and precautions: Wash with water if stabilizing liquid comes in contact with eyes or skin. Do NOT ingest. See MSDS at www.dnagenotek.com.

Small cap, choking hazard.

Storage: 15°C / 30°C (59°F / 86°F)

Summary and explanation of the kit:

OMNigene™ SALIVA DNA and RNA is a collection kit that provides the materials and instructions for collecting and stabilizing human and microbial DNA and RNA from saliva samples.

Collection precautions:

Do NOT eat, drink, smoke, or chew gum for 30 minutes before giving your sample.

Do NOT remove the plastic film from the funnel lid.

Label legend:


- Consult package insert
- Collect sample by (Use by)
- Catalog number
- Caution, consult instructions for use
- Manufacturer
- Lot number
- 15°C / 30°C Storage instructions
- 59°F / 86°F

USER INSTRUCTIONS


Read all instructions prior to collection.

Procedure:


Most people take between 2 and 5 minutes to deliver a sample following steps 1 to 5.

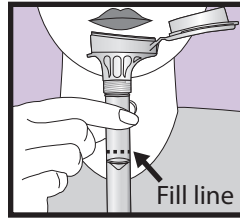


1 Spit into funnel until the amount of liquid (not bubbles) reaches the fill line shown in picture #1.

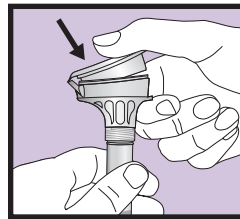


2 Hold the tube upright with one hand. Close the funnel lid with the other hand (as shown) by firmly pushing the lid until you hear a loud click. The liquid in the lid will be released into the tube to mix with the sample. Make sure that the lid is closed tightly.

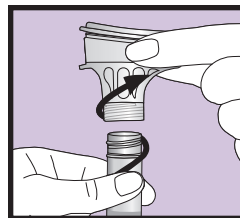

3 Hold the tube upright. Unscrew the funnel from the tube.

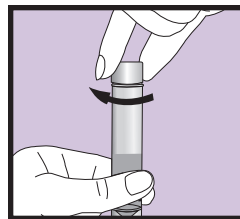

4 Use the small cap to close the tube tightly.

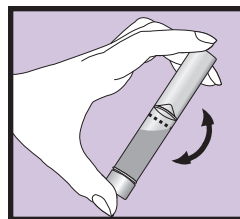

5 Shake the capped tube for 5 seconds. Discard or recycle the funnel.

Collected specimen is potentially infectious and should be handled with appropriate biosafety practices.

Ship in accordance with applicable regulations covering transport of biological specimens.

For Research Use Only
Not for use in diagnostic procedures

DNAGENOTEK™

*Superior samples
Proven performance*

Made in Canada
DNA Genotek Inc.
3000 - 500 Palladium Drive
Ottawa, ON, Canada K2V 1C2

Toll-free (North America): 1.866.813.6354
Tel: +1.613.723.5757 • Fax: +1.613.723.5057
info@dnagenotek.com
www.dnagenotek.com

OMNigene-SALIVA DNA and RNA (OMR-610) is for research use only, not for use in diagnostic procedures. OMNigene and DNA Genotek are trademarks of DNA Genotek, Inc. Some DNA Genotek products may not be available in all geographic regions; contact your sales representative for details. All DNA Genotek protocols, white papers and application notes are available in the support section of our website at www.dnagenotek.com.

Patent (www.dnagenotek.com/legalnotices)
© 2023 DNA Genotek Inc., a subsidiary of OraSure Technologies, Inc., all rights reserved.
PD-PR-01362 Issue 1/2023-02