Liquefaction Reagent protocol

DNAGENOTEK™

www.dnagenotek.com

Tel.: +1.613.723.5757
support@dnagenotek.com
sales@dnagenotek.com

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OMNIgene™ Liquefaction Reagent protocol

Intended use/purpose

OMNIgene™ Liquefaction Reagent (OM-LQR) is intended to liquefy, or render pipettable, stool samples collected in OMNIgene™•GUT that are not pipettable manually and/or by liquid handlers. It is a lab-added reagent used during sample processing to render samples pipettable.

In-use stability statement

Discard the reagent after 12 months of opening or when shelf life is reached — whichever comes first.

Features

- Optimized chemistry for maximum recovery of DNA from stool samples collected with OMNIgene™•GUT.
- Proven to provide consistent results with high molecular weight DNA.
- Scalable purification method for large or small sample volumes.
- Convenient workflow with complete technical support from collection through extraction.
- Cost-effective method that requires minimal equipment.

Materials

- OM-LQR-400 (200 mL) and/or OM-LQR-1600 (800 mL)
- OMNIgene™ Liquefaction Reagent protocol

Warnings and precautions

- For Laboratory Use only.
- Do NOT ingest liquid reagent.
- Do NOT use if packaging is damaged or seal in the funnel lid/cap is broken or leaking.
- Do NOT use OMNIgene™ Liquefaction Reagent beyond the “Use by” date indicated on the reagent bottle.
- Wash with water if reagent comes in contact with eyes or skin. Do NOT ingest.
- Report any serious incident to DNA Genotek and the competent authority in your country.
- Ensure the bottle is tightly capped after each use.
- Refer to SDS for safe disposal of unused reagent.
- Safety Data Sheet (SDS) is available at www.dnagenotek.com.

Product use limitations

Use OMNIgene™ Liquefaction Reagent only as directed in this protocol.

Transportation of OMNIgene™ Liquefaction Reagent

OMNIgene™ Liquefaction Reagent can be transported at room temperature as a laboratory reagent. No special handling is required.

Storage of OMNIgene™ Liquefaction Reagent (Shelf life)

Store at room temperature. Shelf life for OM-LQR-400 (200 mL) and OM-LQR-1600 (800 mL) shall be 30 months when properly capped and stored at room temperature.

Disposal

Discard unused, damaged or leaking kits in accordance with appropriate local, state and federal regulations. Discard as laboratory waste.

Maintenance/repairs

Not applicable. OMNIgene™ Liquefaction Reagent is a reagent — no maintenance or repair required.

Summary of performance characteristics

OMNIgene™ Liquefaction Reagent-purified genomic DNA from an OMNIgene™•GUT stool collection kit provides high-quality and high-quantity DNA sufficient for use in downstream applications, such as PCR, microarray and next generation sequencing.

Product presentations

OMNIgene™ Liquefaction Reagent is available in 2 formats (200 mL and 800 mL) to accommodate the number of preparations required. For example:

Product reference/ Catalog number	Volume of reagent required per OMNIgene™•GUT sample	Number of OMNIgene™•GUT samples that can be prepared
OM-LQR-400 (200 mL bottle)	500 µL	400 samples
OM-LQR-1600 (800 mL bottle)		1600 samples

Warranties

Full terms and conditions for all DNA Genotek products are at <http://www.dnagenotek.com/ROW/terms/index.html>.

Troubleshooting

Contact DNA Genotek technical support at support@dnagenotek.com or call +1 (613) 723-5757, option 6.

Supplied reagent

OMNIgene™ Liquefaction Reagent (DNA Genotek™, Cat. No. OM-LQR-400, OM-LQR-1600)

Note: Perform a visual inspection to assess if precipitation has occurred during storage/shipping; holding bottle against a light is recommended.

If precipitate is present:

1. Incubate at 50°C for 60 minutes.
2. Vigorously shake the bottle for 1 minute by hand. Perform a visual inspection.

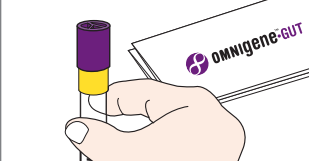
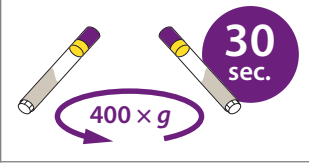
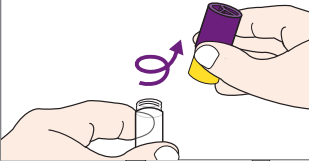
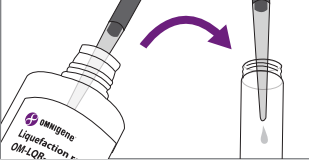
• **Small bottle (OM-LQR-400):** If precipitate is still present, perform additional 50°C incubation for 60 minutes, followed by vigorous shaking for 1 minute by hand. Repeat as necessary until no precipitate remains.

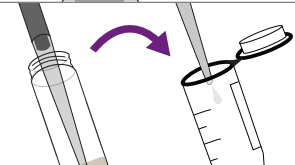
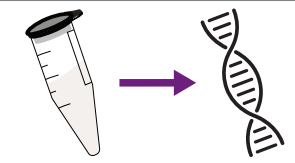
- **Large bottle (OM-LQR-1600):** If precipitate is still present, add a clean 2-inch magnetic stir bar (e.g., VWR®, Cat. No. 58949-038) into the bottle and agitate on a stir plate at 60 rpm for 10 minutes at room temperature. Remove magnetic stir bar after mixing.

Equipment supplied by user

- Pipettors and wide-bore 1000 µL pipette tips (e.g., Axygen™, Cat. No. TF-1005-WB-R-S)
- Bench top vortex mixer
- Bench top centrifuge capable of 400 × g with 15 mL Falcon™ tube -compatible buckets

Procedure

Sample preparation steps	
	1. Collect your fecal sample according to the OMNIgene™•GUT instructions (OM-200: see PD-PR-00612; OMR-200: see PD-PR-01280).
	2. Vortex the sample for 20 seconds.
	3. Centrifuge the sample(s) at 400 × g for 30 seconds. This will collect the material at the bottom of the tube.
	4. With the purple cap still screwed on, unscrew the yellow portion of the tube top and set aside on a clean surface.
	5. Add 500 µL of OMNIgene™ Liquefaction Reagent to the sample. Screw the yellow/purple cap back onto the tube.

Sample preparation steps	
	6. Vortex the sample for 60 seconds. Sample will become visibly homogenous and liquid once the reagent is fully mixed in. IMPORTANT: If you are not proceeding immediately to extraction, repeating step 6 prior to proceeding to step 7 is required to guarantee a homogenous sample.
	7. With the purple cap still screwed on, unscrew the yellow portion of the tube top and set aside on a clean surface. Using a wide-bore 1000 µL pipette tip, slowly pipette the sample into the extraction tube/vessel. IMPORTANT: Slow aspiration is critical.
	8. Proceed with preferred extraction method.

Label legend:

IVD

In vitro diagnostic medical device

UK

UKCA marking

REF

Catalog number

CE

CE marking

Manufacturer

Consult electronic instructions for use

Caution

EC REP

European Authorized Representative

CH REP

Swiss Authorized Representative

LOT

Lot number

UDI

Unique Device Identifier

QTY

Quantity of units per package

Use by/In-use stability

15°C / 25°C
59°F / 77°F

Storage instructions

MADE IN US
FROM FOREIGN
AND DOMESTIC
MATERIALS

Made in US from foreign
and domestic materials

Technical support is available Monday to Friday (9h00 to 17h00 ET):

Toll-free (North America): 1.866.813.6354, option 6
All other countries: +1.613.723.5757, option 6
Email: support@dnagenotek.com

DNA Genotek Inc.
3000 - 500 Palladium Drive
Ottawa, ON, K2V 1C2, Canada

UK Responsible Person: Qarad UK Ltd., 8 Northumberland Ave, Westminster, London, WC2N 5BY United Kingdom
EC REP Qarad EC-REP BV, Pas 257, 2440 Geel, Belgium
CH REP Qarad Suisse S.A., World Trade Center, Avenue Gratta-Paille 2, 1018 Lausanne, Switzerland
Australian Sponsor: Emergo Australia, Level 20, Tower II, Darling Park, 201 Sussex Street, Sydney, NSW 2000 Australia

Some DNA Genotek products may not be available in all geographic regions; contact your sales representative for details.
OMNigene and DNA Genotek are trademarks of DNA Genotek Inc., registered under various jurisdictions.
All other brands and names contained herein are the property of their respective owners.
All DNA Genotek protocols, white papers and application notes are available in the support section of our website at www.dnagenotek.com.
Patent (www.dnagenotek.com/legalnotices)

© 2024 DNA Genotek Inc., a subsidiary of OraSure Technologies, Inc., all rights reserved.

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Odgovorna oseba v ZK: Qarad UK Ltd., 8 Northumberland Ave, Westminster, London, WC2N 5BY Združeno kraljestvo
CH REP Qarad Suisse S.A., World Trade Center, Avenue Gratta-Paille 2, 1018 Lausanne, Švica
OMNIGENE in DNA Genotek sta blagovni znamki DNA Genotek Inc.
© 2024 DNA Genotek Inc., hčerinska družba OraSure Technologies, Inc., vse pravice pridržane.