
For Research Use Only
Not for use in diagnostic procedures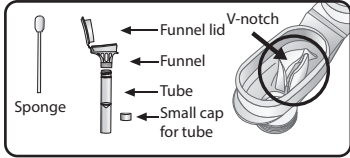

Read all instructions prior to collection.

Intended use: This product is designed for the assisted collection of human DNA from saliva samples.

Contents: 1 kit containing 0.75 mL stabilizing liquid

Warnings and precautions:

- Choking hazards:
 - Small cap in collection kit.
 - Plastic bag containing sponges.
 - Caution should be used when inserting sponge into donor's mouth.
- For supervised collections:
 - Do NOT leave donor unattended.
 - Do NOT allow donor to handle the sponge, small cap or packaging.
 - Wash with water if stabilizing liquid comes in contact with eyes or skin. Do not ingest. See SDS at www.dnagenotek.com.

Storage: 15°C / 30°C (59°F / 86°F)

Summary and explanation of the kit:

Oragene™-DISCOVER is an assisted collection kit that provides the materials and instructions for collecting and stabilizing saliva specimens.

Collection precautions:

Do NOT remove the plastic film from the funnel lid.

Check sponge for damage each time before inserting into donor's mouth. Use second sponge if first sponge shows any signs of wear or tear.

Do NOT substitute with other sponges or swabs.

Procedure:

- Ensure donor does NOT eat, drink, smoke, or chew gum for 30 minutes before collecting a saliva sample.
- Ensure donor is in an upright position during sample collection.
- It may take up to 15 minutes to collect a saliva sample following steps 1 to 7.

DNagenotek™

*Superior samples
Proven performance*

■ DNA Genotek Inc.
3000 - 500 Palladium Drive
Ottawa, ON, Canada
K2V 1C2

Tel.: +1.613.723.5757
Fax: +1.613.723.5057
info@dnagenotek.com
www.dnagenotek.com


PD-PR-619 v2


ENGLISH

Place one sponge in cheek pouch. Gently move the sponge along the gums and inner cheeks for 30 seconds to soak up as much saliva as possible.

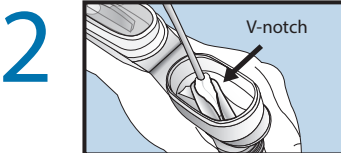
Once sponge is saturated with saliva, insert sponge in V-notch of funnel. Wring saliva out of sponge using a twisting and pushing motion against the inner wall of the V-notch. Saliva will flow into tube.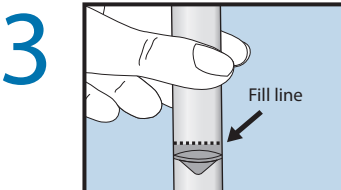


Repeat these steps (1 to 2) USING THE SAME SPONGE until the liquid saliva (not bubbles) reaches the fill line. Check sponge for damage each time before inserting into donor's mouth. Use second sponge if first sponge shows any signs of wear or tear.


Tap tube bottom against hard surface to reduce number of bubbles.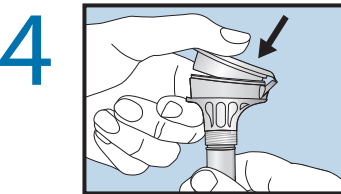


Hold the tube upright with one hand. Close the lid with the other hand (as shown) by firmly pushing the lid until you hear a loud click. The liquid in the lid will be released into the tube to mix with the saliva. Make sure the lid is closed tightly.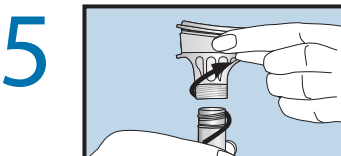

Hold the tube upright. Unscrew the funnel from the tube.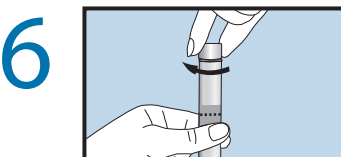

Use the small cap to close the tube tightly.

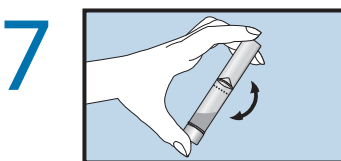Shake the capped tube for 5 seconds. Discard or recycle the funnel. Discard sponges.

REF OGR-575 and REF OGR-675

Oragene is a trademark of DNA Genotek Inc., registered in various jurisdictions. Some DNA Genotek products may not be available in all geographic regions; contact your sales representative for details. All DNA Genotek protocols, white papers and application notes are available in the support section of our website at www.dnagenotek.com.

Patent (www.dnagenotek.com/legal-notice)
© 2025 DNA Genotek Inc., a subsidiary of OraSure Technologies, Inc., all rights reserved.

Label legend:

-  Flammable liquid
-  Consult package insert
-  Collect saliva by (Use by)
-  Catalog number
-  Caution, consult instructions for use
-  Storage instructions
-  Manufacturer
-  Do not re-use
-  Lot number

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© 2025 DNA Genotek Inc., ein Tochterunternehmen von OraSure Technologies, Inc., alle Rechte vorbehalten.