


For Research Use Only
Not for use in diagnostic procedures

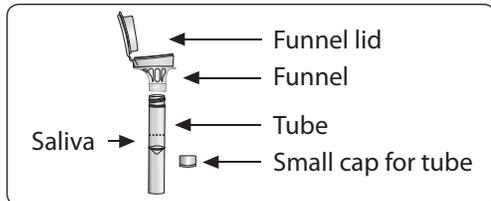

Collection precautions:

Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.

Do NOT remove the plastic film from the funnel lid.

This product is designed for the collection of human DNA from saliva samples.

Contents: Kit contains stabilizing liquid.

Warnings and precautions: Wash with water if stabilizing liquid comes in contact with eyes or skin. Do not ingest. See MSDS at www.dnagenetek.com.

Small cap, choking hazard.

Storage: 15°C / 30°C

Summary and explanation of the kit:

Oragene-DISCOVER is a self-collection kit that provides the materials and instructions for collecting and stabilizing saliva specimens.

U.S. Patent Nos. 8,221,381 and 7,482,116;
European Patent No. 1 513 952 and 1 956 969; Patent pending
Canadian Design Nos. 127470; 132896; 132897
U.S. D631,554 S and D640,795 S
Community Design Nos. 001095186-0001; -0002; -0003

*Oragene is a registered trademark of DNA Genetek Inc.
Oragene®-DISCOVER is for research use only, not for use in diagnostic procedures.
Some DNA Genetek products may not be available in all geographic regions.


PD-PR-00279 Issue 1/2012-10
© 2012 DNA Genetek Inc., a subsidiary of OraSure Technologies, Inc., all rights reserved.

USER INSTRUCTIONS


Read all instructions prior to collection

Procedure:


Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 5.

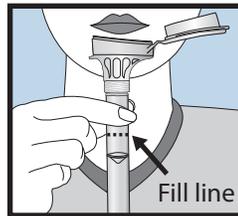


- 1 Spit into funnel until the amount of liquid saliva (not bubbles) reaches the fill line shown in picture #1.

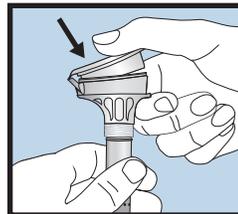


- 2 Hold the tube upright with one hand. Close the funnel lid with the other hand (as shown) by firmly pushing the lid until you hear a loud click. The liquid in the lid will be released into the tube to mix with the saliva. Make sure that the lid is closed tightly.

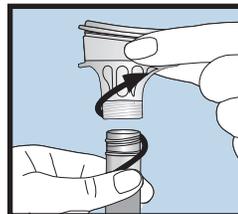

- 3 Hold the tube upright. Unscrew the funnel from the tube.

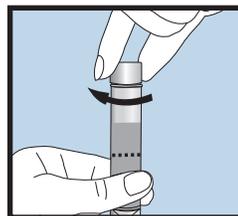

- 4 Use the small cap to close the tube tightly.

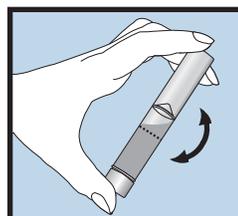

- 5 Shake the capped tube for 5 seconds. Discard or recycle the funnel.

Label legend:

	Consult package insert
	Collect saliva by (Use by)
	Catalog number
	Industrial Design Patent
	Caution, consult instructions for use
	Storage instructions
	Manufacturer

For Research Use Only
Not for use in diagnostic procedures

Superior samples • Proven performance

Made in Canada by DNA Genetek Inc.
2 Beaverbrook Road, Ottawa, ON, Canada K2K 1L1
Toll-free (North America): 1.866.813.6354
Tel: 613.723.5757 • Fax: 613.723.5057
www.dnagenetek.com • info@dnagenetek.com

DNAgenetek



www.dnagenetek.com