


Oral sample collection instructions

- Animal should NOT eat for 30 minutes or drink for 10 minutes before saliva collection.
- Do NOT scrape the animal's teeth or cheek with the sponge.
- Do NOT allow the animal to chew or bite the sponge.

Precautions:


Tube contains stabilizing liquid. Do NOT spill. This liquid preserves the DNA sample and is required by the lab to analyze the sample.

IMPORTANT: Do NOT remove the cap prior to sample collection.

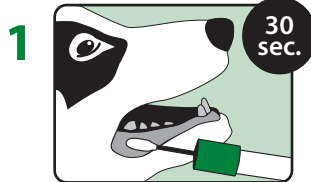

- Place the sponge in animal's mouth at the cheek pouch as shown. Collect saliva for **30 seconds** by moving sponge and mopping saliva where it naturally pools (in the cheek pouch and under the tongue).
 - For animals older than 6 months, moderate restraint may be required.


IMPORTANT: Tube contains stabilizing liquid. Do NOT spill.

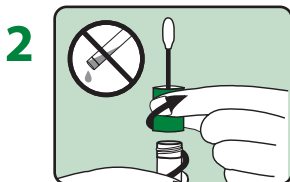


- Hold the tube upright and unscrew the cap from the tube.



- Turn the cap upside down and place the sponge in the tube.
 - Screw cap on tightly to prevent sample from leaking during transport.

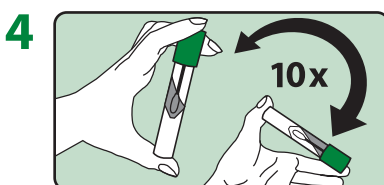

- Invert and shake the tube vigorously **10 times** to thoroughly mix sample.

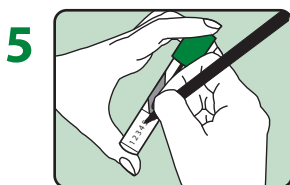

- Optional: Using a permanent marker, clearly write the animal identification number on the white space available on the tube label.

Intended use: This product is designed for the collection of DNA from animals.

Warnings: Wash with water if stabilizing liquid comes in contact with eyes or skin. Do NOT ingest.

Storage: 15°C-25°C (59°F-77°F)

Material Safety Data Sheet (MSDS) available at www.dnagenotek.com

For Animal Use Only

Patent (www.dnagenotek.com/legalnotices)
 PD-PR-00347 Issue 3/2021-09
 © 2021 DNA Genotek Inc., all rights reserved.

DNA Genotek Inc.
 3000 - 500 Palladium Drive
 Ottawa, ON, Canada K2V 1C2
 Subsidiary of OraSure Technologies, Inc.

Toll-free (North America): 1.866.813.6354
 Tel.: 613.723.5757 • Fax: 613.723.5057
www.dnagenotek.com
info@dnagenotek.com

DNAGENOTEK®

www.dnagenotek.com