



for pediatrics

REF OC-175



Collection precautions:

Ensure the sponge tip does NOT come into contact with any surface prior to collection.

Donor should NOT eat, drink, smoke or chew gum for 30 minutes before collecting oral sample.

Intended for the collection of human DNA from pediatric oral samples.

Contents: Contains 1 collection kit.

Warning and precautions: Sponge tip is a choking hazard. Caution should be used when inserting sponge into the mouth. For assisted collections do NOT leave infant or child unattended.

Wash with water if liquid comes in contact with eyes or skin. Do NOT ingest. See MSDS at www.dnagenotek.com

Storage: 15-25°C

Summary and explanation of the kit:

ORACollect for pediatrics is an assisted collection kit that provides the materials and instructions for collecting and stabilizing human DNA from pediatric oral samples.

Label legend:

- Manufacturer
REF Catalog number
Non-sterile
LOT Lot number
Collect sample by (Use by)
Storage instructions 15°C-25°C

Instructions for sample collection:

5-step instructional diagram for sample collection: 1. Insert sponge into mouth and rub 10x. 2. Repeat on opposite side 10x. 3. Unscrew cap. 4. Insert sponge into cap. 5. Shake tube 10x.

For Research Use Only
Not for use in diagnostic procedures



Superior samples
Proven performance

DNA Genotek Inc.
3000 - 500 Palladium Drive
Ottawa, ON, Canada K2V 1C2

Toll-free (North America): 1.866.813.6354
Tel: +1.613.723.5757
Fax: +1.613.723.5057
info@dnagenotek.com
www.dnagenotek.com

*ORACollect is a registered trademark of DNA Genotek Inc.
Some DNA Genotek products may not be available in all geographic regions, contact your sales representative for details.
All DNA Genotek protocols, white papers and application notes, are available in the support section of our website at www.dnagenotek.com.
Patent (www.dnagenotek.com/legalnotices)
© 2019 DNA Genotek Inc., a subsidiary of OraSure Technologies, Inc., all rights reserved.
PD-PR-00542 Issue 2/2019-10